


# Operating Procedure – MetaVis Microscope

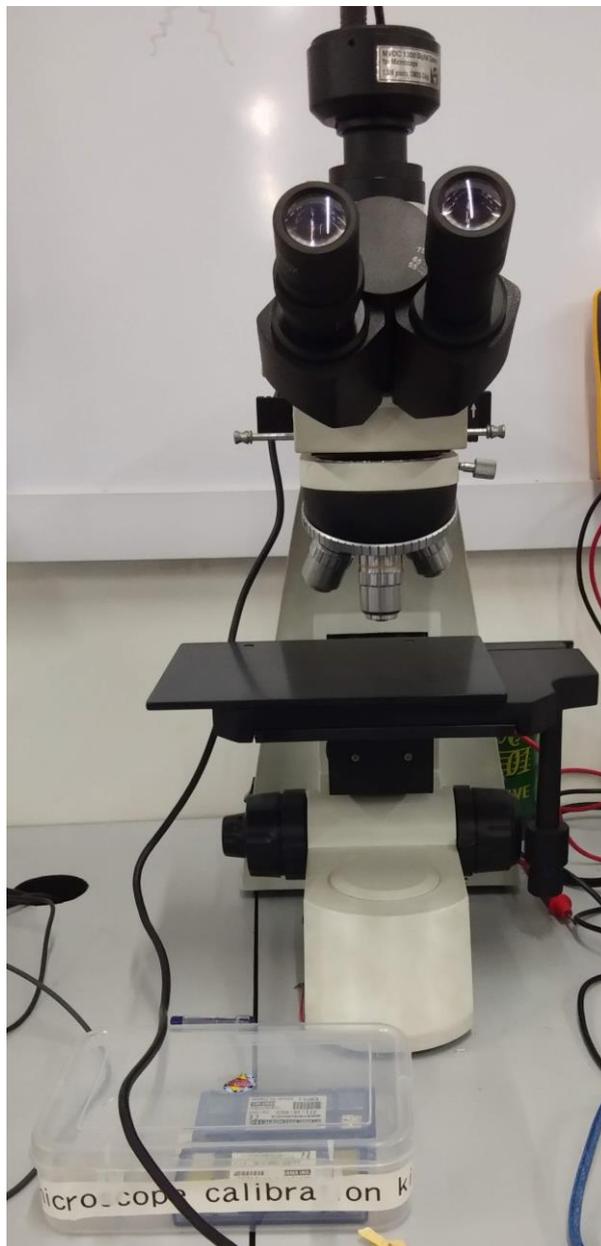

**National Centre for Photovoltaic Research and Education (NCPRE)**

**Indian Institute of Technology- Bombay**

**Powai, Mumbai-400076**

**INDIA**

## **Standard Operating Procedure**

1. Switch on the power supply button on the microscope.
2. Switch on the computer.
3. Open the microscope software (ScopePhoto( $\times 86$ ))
4. Choose proper objective and adjust the desired area and focus by moving the stage in X, Y and Z directions.
5. Capture the image using software (in ImageJ) and save in appropriate folder.
6. Switch off the computer.
7. Switch off the power supply button on the microscope.



**End of Document**